## Change Password in Freezerpro

When you logged in into Freezerpro you can change your password

Home		
	🖂 👻 🛄 👻 Actions 👻 Search: Live	e search re
Bookmarks	About Add New Freezer	
	G Define New Sample Ty	ре
My Samples     My Boxes	<ul> <li>About Fr</li> <li>Terms of G Import Samples</li> </ul>	
User Preferences	<ul> <li>Privacy P G Update</li> </ul>	•
	Support O My Samples	
	<ul> <li>User Guic</li> </ul>	
	FreezerPt Q Advanced Search	
	Ask a Qu     Q     Batch Search     Our Prod	•
	Alerts	
	III Freezers Sensor Monito	or
	Reports And Audit	•
	Browse Samples	•
	8 Browse Sample Groups	5
	Source Browse Sample Source	s
	a User Preferences	

At the Home tab, click on Action menu and choose User Preferences. A pop-up window appears

User Preferences			×	<
Country Settings:	Sweden		•	
Page Size (Samples per page):	100			
Barcode Printer:	ZZ-SMB printer Box Label		•	
After Login Open:	O My Samples List			
	O My Last Viewed Box			
	O Samples Added This Week			
	O No Action			
	Change Password			
		Save	Cancel	

## Click on the Change Password button. A new pop-up window appears

Change Password		×
Old Password: Password: Verify Password:		
	Change	Cancel

Type your password and then type your new password twice. Click on Change button to change the password.